

GlitterBuddy™ Kit



GlitterBuddy™ Kit provides a convenient, fun & effective handwashing and cross contamination training session.

Rating: Not Rated Yet

Price:

Variant price modifier:

Base price with tax:

Price with discount:

Salesprice with discount:

Sales price:

Sales price without tax:

Discount:

Tax amount:

[Ask a question about this product](#)

Description

Made of durable paper-board, the GlitterBuddy™ Kit folds down to a compact box that contains all you need to run a succesful hygiene traininig session for both hand washing and cross contamination.

When unfolded, the GlitterBuddy™ Kit resembles a germ, and you simply look through his mouth to see the results of the training session.

To demonstrate hand washing technique:-

1. Simply rub the GlitterBug® Potion onto your hands like you would with hand cream.
2. Wash your hands as you would normally and dry them.
3. Place your hands inside the GlitterBuddy™ Kit and look thru his mouth under the UV light for any fluorescent areas. These fluorescent areas highlight spots that were not adequately washed during the handwashing procedur - thus areas where microorganisms may like to lurk on your hands.

To demonstrate cross contamination using GlitterBug® Powder :-

1. Sprinkle some powder onto a piece of paper.
2. Pretend to sneeze on the paper.
3. Ask people to take the paper and sign their name for attendance records.
4. Using the GlitterBuddy™ Kit look at the fingers of the first person taking the paper. The glowing area demonstrates the cross contamination of 'bugs' from the paper to their hands.

Training : GlitterBuddy™ Kit

The kit is supplied with:-

- SpotShooter® 8 UV Lamp
- GlitterBug® Potion (50mL)
- GlitterBug® Powder (18g)
- GlitterBug® Instruction Manual
- Hand wash DVD
- How to wash your hands poster
- Two pages of 'I was rude' stickers
- GlitterBug® Magnet
- Two GlitterBug® Hand Brushes



Ordering Details

GlitterBuddy™ Kit

Product Code: BRBUDDYKIT