

Hydrion® Quat Test Papers



Hydrion® Quat Test Papers are an easy, reliable and affordable method for the analysis of quaternary ammonium-based sanitisers in solutions.

Rating: Not Rated Yet

Price

[Ask a question about this product](#)

Description

Hydrion® Sanitiser test papers are used to determine the levels of quaternary ammonium-based sanitiser in solutions. Three products are available:- QT10 is formulated for 2-chain quat compounds with measurements in the range of 0-400ppm; QT-40 is for 4-chain quats with measurements of 0-500ppm; and QC1001 is for high range 2-chain quats with a measurement range of 0-1000ppm.

Instructions:

1. Dip the strip into the sanitising solution for 10 seconds.
2. Instantly compare the resulting colour with the enclosed colour chart.
3. Please note that the test solution should be between 18-25°C.

Storage Conditions:

Room Temperature

Ordering Details:

- | | |
|----------------------|-----------------------|
| ▪ Range: 0 - 400ppm | Product Code: QT-10 |
| ▪ Range: 0 - 500ppm | Product Code: QT-40 |
| ▪ Range: 0 - 1000ppm | Product Code: QC-1001 |

